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Ziegelei 1

D-72336 Balingen

E-Mail: info@kern-sohn.com

Tel: +49-[0]7433- 9933-0

Fax: +49-[0]7433-9933-149

Internet: www.kern-sohn.com

Betriebs- und Installationsanleitung Anzeigegeräte
Operating and Installation Instructions Display units
Notice d'utilisation et d'installation Afficheurs

KERN VHP-T

Version 1.1

05/2016

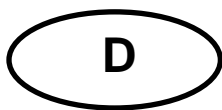
Deutsch

English

Français

VHP-T-BA_IA-def-1611

- D** Weitere Sprachversionen finden Sie online unter www.kern-sohn.com/manuals
- CZ** Další jazykové verze najdete na webu pod adresou www.kern-sohn.com/manuals
- DK** Yderligere sprogversioner finder de online på www.kern-sohn.com/manuals
- E** Más versiones de idiomas se encuentran online bajo www.kern-sohn.com/manuals
- EST** Rohkem keeli internetis aadressil www.kern-sohn.com/manuals
- F** Vous trouverez d'autres versions de langue online sous www.kern-sohn.com/manuals
- GB** Further language versions you will find online under www.kern-sohn.com/manuals
- H** A használati utasítás egyéb nyelveken a www.kern-sohn.com/manuals címről tölthető le
- I** Trovate altre versioni di lingue online in www.kern-sohn.com/manuals
- N** Ytterligere språkversjoner finner du online under www.kern-sohn.com/manuals
- NL** Bijkomende taalversies vindt u online op www.kern-sohn.com/manuals
- P** Encontram-se online mais versões de línguas em www.kern-sohn.com/manuals
- PL** Inne wersje językowe znajdują Państwo na stronie www.kern-sohn.com/manuals
- S** Ytterligare språkversioner finns online under www.kern-sohn.com/manuals
- SF** Muita kieliversioita löydät osoitteesta www.kern-sohn.com/manuals
- SLO** Ostale jezikovne različice boste našli online na www.kern-sohn.com/manuals
- TR** Diğer lisan versiyonlarını internetten www.kern-sohn.com/manuals adresinden temin edebilirsiniz

KERN VHP-T

Version 1.1 05/2016

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KERN VHP-T

Version 1.1 05/2016

Operating and installation instructions Display units

Contents

1	Technical data	4
2	Device overview VHP-T	5
3	Keyboard overview	6
3.1	Overview of display	7
4	Basic Information (General)	7
4.1	Proper use	7
4.2	Improper Use	7
4.3	Warranty	8
4.4	Monitoring of Test Resources	8
5	Basic Safety Precautions	8
5.1	Pay attention to the instructions in the Operation Manual.....	8
5.2	Personnel training.....	8
6	Transport and storage	9
6.1	Testing upon acceptance	9
6.2	Packaging / return transport	9
7	Unpacking and placing	10
7.1	Installation Site, Location of Use	10
7.2	Unpacking.....	11
7.3	Scope of delivery / serial accessories:	11
7.4	Placing	11
7.5	Storage battery operation (optional).....	11
7.6	Selection of platform quantity	12
7.7	Adjustment.....	12
7.8	How to carry out adjustments	14
8	Operation	15
8.1	Start-up	15
8.2	Switching Off	15
8.3	Zeroing	15
8.4	Simple weighing	15
8.5	Weighing with tare	16
8.6	Manual totalizing.....	17
8.7	Display background illumination.....	17
8.8	Automatic switch-off function „AUTO OFF“	18

9	Menu	19
10	Servicing, maintenance, disposal	20
10.1	Cleaning	20
10.2	Servicing, maintenance	20
10.3	Disposal	20
10.4	Error messages	20
11	Instant help.....	21
12	Configure display unit	22
13	Declaration of Conformity.....	23

1 Technical data

KERN	VHP-T	
Display	5-digit	
Resolution	3000	
Divisions	1,2,5,10,20,50	
Weighing Units	Kg	
Functions	Totalization	
Display	LCD 10 mm digits with back lighting	
DMS weighing cells	Max. 6 items 700 Ω; sensitivity: 1±0.2mv/V.	
Range calibration	We recommend ≥ 50 % max.	
Data output	RS232	
Electric Supply	Input voltage 100 V – 240 V	
	Installed rechargeable battery: DC 6V 10Ah	
Housing	With case: 510x340x220	Without case: 350x280x170
Admissible ambient temperature	0°C – 40°C	
Net weight	6 kg	
Operating / charge time of battery	13 h / 12 h	
RS 232 interface	Serially	
IP protection	IP 54	

2 Device overview VHP-T

1. Printer
2. Display
3. Keyboard
4. Connections (platforms, load connection of battery)
5. Case

3 Keyboard overview

Button	Function
	<ul style="list-style-type: none"> • Switch-on
	<ul style="list-style-type: none"> • Switch-off
	<ul style="list-style-type: none"> • Taring
	<ul style="list-style-type: none"> • Display sum total
	<ul style="list-style-type: none"> • Store weighing value
	<ul style="list-style-type: none"> • Delete weighing values
        	<ul style="list-style-type: none"> • Printing • Return key • Setting the date • Setting the time • Entry of license plate • Switch on and off background illumination • Input of letters • Input of numbers • Zero setting key

3.1 Overview of display

Display	Significance
	Battery very low
G	Gross weight
NT	Net weight
Kg	Weighing unit

4 Basic Information (General)

4.1 Proper use

The display unit acquired by you is used in combination with a weighing plate and serves to determine the weighing value of material to be weighed. It is intended to be used as a “non-automatic weighing system”, i.e. the material to be weighed is manually and carefully placed in the centre of the weighing plate. As soon as a stable weighing value is reached the weighing value can be read.

4.2 Improper Use

Do not use display unit for dynamic weighing. In the event that small quantities are removed or added to the material to be weighed, incorrect weighing results can be displayed due to the “stability compensation” in the display unit. (Example: Slowly draining fluids from a container on the balance.)

Do not leave permanent load on the weighing pan. This may damage the measuring system.

Impacts and overloading exceeding the stated maximum load (max) of the weighing plate, minus a possibly existing tare load, must be strictly avoided. Both, the weighing plate and the display unit may be damaged during this process.

Never operate display unit in explosive environment. The serial version is not explosion protected.

Changes to the display unit's design are not permitted. This may lead to incorrect weighing results, safety-related faults and destruction of the display unit.

The display unit may only be operated in accordance with the described default settings. Other areas of use must be released by KERN in writing.

4.3 Warranty

Warranty claims shall be voided in case

- Our conditions in the operation manual are ignored
- The appliance is used outside the described uses
- The appliance is modified or opened
- Mechanical damage or damage by media, liquids, natural wear and tear
- The appliance is improperly set up or incorrectly electrically connected
- The measuring system is overloaded

4.4 Monitoring of Test Resources

In the framework of quality assurance the measuring-related properties of the display unit and, if applicable, the testing weight, must be checked regularly. The responsible user must define a suitable interval as well as type and scope of this test. Information is available on KERN's home page (www.kern-sohn.com) with regard to the monitoring of display units' test substances and the test weights required for this. In KERN's accredited DKD calibration laboratory test weights and display units may be calibrated (return to the national standard) fast and at moderate cost.

5 Basic Safety Precautions

5.1 Pay attention to the instructions in the Operation Manual

- ⇒ Carefully read this operation manual before setup and commissioning, even if you are already familiar with KERN balances.
- ⇒ All language versions contain a non-binding translation. The original German is binding.

5.2 Personnel training

The appliance may only be operated and maintained by trained personnel.

6 Transport and storage

6.1 Testing upon acceptance

When receiving the appliance, please check packaging immediately, and the appliance itself when unpacking for possible visible damage.

6.2 Packaging / return transport

- ⇒ Keep all parts of the original packaging for a possibly required return.
- ⇒ Only use original packaging for returning.
- ⇒ Prior to dispatch disconnect all cables and remove loose/mobile parts.
- ⇒ Reattach possibly supplied transport securing devices.
- ⇒ Secure all parts such as the glass wind screen, the weighing platform, power unit etc. against shifting and damage.

7 Unpacking and placing

7.1 Installation Site, Location of Use

The display units are designed in a way that reliable weighing results are achieved in common conditions of use.

Precise and fast work is achieved by selecting the right place for your display unit and your weighing plate.

On the installation site observe the following:

- Place the display unit and the weighing plate on a stable, even surface.
- Avoid extreme heat as well as temperature fluctuation caused by installing next to a radiator or in the direct sunlight;
- Protect the display unit and the weighing plate against direct draft from open windows or doors.
- Avoid jarring during weighing;
- Protect the display unit and the weighing plate against high humidity, vapours and dust.
- Do not expose the display unit to extreme dampness for longer periods of time. Non-permitted condensation (condensation of air humidity on the appliance) may occur if a cold appliance is taken to a considerably warmer environment. In this case, acclimatize the disconnected appliance for ca. 2 hours at room temperature.
- Avoid static charge of goods to be weighed or weighing container.

Major display deviations (incorrect weighing results) may be experienced should electromagnetic fields (e.g. due to mobile phones or radio equipment), static electricity accumulations or instable power supply occur. Change location or remove source of interference.

7.2 Unpacking

Carefully remove the display unit from packaging, remove plastic cover and place it in the designated work area.

and place it in the designated work area.

7.3 Scope of delivery / serial accessories:

- Display Unit
- Mains adapter for loading the battery
- RS 232 interface cable
- Operating manual
- 1 x paper roll for printer

7.4 Placing

Mount the display unit in a way that facilitates operation and where it is easy to see.

7.5 Storage battery operation (optional)

Before the first use, the battery should be charged by connecting it to the mains power supply for at least 12 hours.

Charge the battery with the help of the supplied power pack.

The LED display at the mains adapter informs you during loading about the loading status of the rechargeable battery.

red: Voltage has dropped below prescribed minimum.

green: Rechargeable battery completely reloaded

To conserve energy, enable the automatic switch-off function „AUTO OFF“, see chap. 8.8.

7.6 Selection of platform quantity

1.	Turn on device using the  button
2.	Use  button to change into the window for date input.
3.	Now enter the following password: 878586 (the last two ciphers can be ignored)
4.	Use the C button to change between the individual parameters.
5.	Select the parameter (8): Mode
6.	Confirm by pressing the  button
7.	Now enter the quantity of platforms. As per standard there are 2. (up to 6 platforms available)

7.7 Adjustment

As the acceleration value due to gravity is not the same at every location on earth, each display unit with connected weighing plate must be coordinated - in compliance with the underlying physical weighing principle - to the existing acceleration due to gravity at its place of location (only if the weighing system has not already been adjusted to the location in the factory). This adjustment process must be carried out for the first commissioning, after each change of location as well as in case of fluctuating environment temperature. To receive accurate measuring values it is also recommended to adjust the display unit periodically in weighing operation.

i	<ul style="list-style-type: none">• Prepare the required adjustment weight. The weight to be used depends on the capacity of the scale. Carry out adjustment as near as possible to the scale's maximum weight. Info about test weights can be found on the Internet at: http://www.kern-sohn.com.• Observe stable environmental conditions. Stabilisation requires a certain warm-up time.
----------	---

Call up menu:

1.	Turn on device using the  - button
2.	Use  button to change into the window for date input.
3.	Now enter the following password: 878586 (the last two ciphers can be ignored)
4.	Use the C button to change between the individual parameters.
5.	Select the parameter (5): Select Calibration
6.	Confirm by pressing the  button

7.8 How to carry out adjustments


Important: Always adjust both weighing platforms in a parallel manner. Not one by one. Always place the adjustment weight evenly on both platforms. In menu define the respective platform as a half.

Example: Place 15 tons of adjustment weight on both platforms. In the menu adjust each platform with the following input:

Platform 1: 7500kg

Platform 2: 7500kg

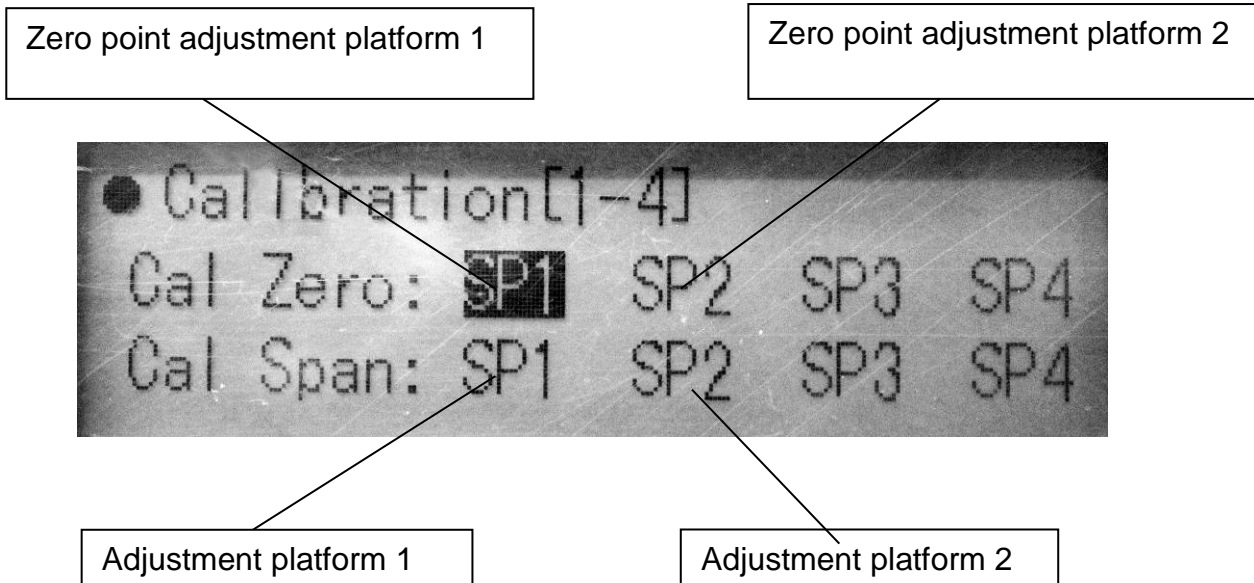The following display appears after acknowledgement of parameter (5) Calibration:

⇒ To select the platforms, press the C button

⇒ Use the  button to select the respective platform.

⇒ Carry out zero point adjustment (CAL ZERO) at platform 1 (SP1) and platform 2 (SP2). During zero point adjustment the weighing plate must not be charged.

⇒ Always place the adjustment weight evenly on both platforms.

⇒ Carry out adjustment at platform 1 (SP1) and platform 2 (SP2). (please note the example above: 50% of the total load per platform must be defined.)

8 Operation

8.1 Start-up

- ⇒ Press . As soon as the weight display appears, the instrument will be ready to weigh.

8.2 Switching Off

- ⇒ Press , the display will disappear.

8.3 Zeroing

Resetting to zero corrects the influence of light soiling on the weighing plate. If required, the device may be reset to zero at any time.

- ⇒ To unload the weighing system

- ⇒ Press , the zero display appears.

8.4 Simple weighing

- ⇒ Place/drive-on weighing goods (at walking speed)
⇒ Read weighing result.

Overload warning

Overloading exceeding the stated maximum load (max) of the device, minus a possibly existing tare load, must be strictly avoided. This could damage the instrument.

Exceeding the maximum load is indicated by the display „----“.

Unload weighing system or reduce preload.

8.5 Weighing with tare

- ⇒ Deposit weighing vessel. After successful standstill control press the

The zero display appears.

The weight of the container is now internally saved.

- ⇒ Weigh the material, the net weight will be indicated.
- ⇒ The weight of the weighing container will be displayed as a minus number after removing the weighing container.
- ⇒ The tare procedure can be repeated as many times as necessary, for example with initial weighing of several components for a mix (add-on weighing). The limit is reached when the taring range reaches its capacity limit.

- ⇒ To delete the tare value, remove load from weighing plate and press

8.6 Manual totalizing

 PRINT

After pressing the  button, the separate weighing results of the platforms are added up to a total result (axle) and printed out.

See printout example below:

Axle load - report	
30/11/2014	16:03:34
SN. 2	
HGV license plate:	
BLM	
Article: 01	

Wheel 1:	70 kg
Wheel 2:	40 kg

Axle 1:	110 kg

G =	110 kg
T =	0 kg
NT=	110 kg
User:	

8.7 Display background illumination

⇒ The background illumination can be turned on and off by using  button.

8.8 Automatic switch-off function „AUTO OFF“

1.	Turn on device using the  - button
2.	Use  button to change into the window for date input.
3.	Now enter the following password: 878586 (the last two ciphers can be ignored)
4.	Use the C button to change between the individual parameters.
5.	Select the parameter (9): Select "Auto-Off time".
6.	Confirm by pressing the  button
7.	Now a time of 1-99 minutes can be defined.
8.	Confirm by pressing the  button

9 Menu

Navigation in the menu:

Access to menu

1.	Turn on device using the  button
2.	Use  button to change into the window for date input.
3.	Now enter the following password: 878586 (the last two ciphers can be ignored)
4.	Use the C button to change between the individual parameters.
5.	Using  button select the corresponding parameter.

Menu overview

Menu	Function
(1) Capacity	<ul style="list-style-type: none">Setting weighing scale capacity
(2) Readability	<ul style="list-style-type: none">Setting weighing scale readability.
(3) Decimal point	<ul style="list-style-type: none">Setting decimal point (decimal place).
(4)	<ul style="list-style-type: none">(please do not modify standard setting on „5“)
(5) Calibration	<ul style="list-style-type: none">Adjustment of balance.
(6)	<ul style="list-style-type: none">Not documented
(7) Language	<ul style="list-style-type: none">Selection of language German and English
(8) Mode	<ul style="list-style-type: none">Selection of quantity of platforms (up to 6 platforms possible)

10 Servicing, maintenance, disposal

10.1 Cleaning

- Before cleaning, disconnect the appliance from the operating voltage.
- Do not use aggressive detergents (solvents or similar).

10.2 Servicing, maintenance

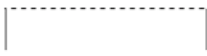
The appliance may only be opened by trained service technicians who are authorized by KERN.

Before opening, disconnect from power supply.

10.3 Disposal

Disposal of packaging and appliance must be carried out by operator according to valid national or regional law of the location where the appliance is used.

10.4 Error messages

Error message	Description	Possible causes
+/-RRR(EEE)	No defined signal	<ul style="list-style-type: none">• Weighing platform not connected.
	Overload display	<ul style="list-style-type: none">• Too much weight on the weighing platform.
	Underload display	<ul style="list-style-type: none">• Insufficient weight on the weighing platform.

Should other error messages occur, switch balance off and then on again. If the error message remains inform manufacturer.

11 Instant help

In case of an error in the program process, briefly turn off the display unit and disconnect from power supply. The weighing process must then be restarted from the beginning.

Help:

Fault

Possible cause

The displayed weight does not glow.

- The display unit is not switched on.
- Mains power supply interrupted (mains cable defective).
- Power supply interrupted.
- (Rechargeable) batteries are inserted incorrectly or empty
- No (rechargeable) batteries inserted.

The displayed weight is permanently changing

- Draught/air movement
- Table/floor vibrations
- Weighing pan has contact with other objects.
- Electromagnetic fields / static charging (choose different location/switch off interfering device if possible)

The weighing result is obviously incorrect

- The display of the balance is not at zero
- Adjustment is no longer correct.
- Great fluctuations in temperature.
- Warm-up time was ignored.
- Electromagnetic fields / static charging (choose different location/switch off interfering device if possible)

Should other error messages occur, switch display unit off and then on again. If the error message remains inform manufacturer.

12 Configure display unit

1.	Turn on device using the  - button
2.	Use  button to change into the window for date input.
3.	Now enter the following password: 878586 (the last two ciphers in the display can be ignored)
4.	Use the C button to change between the individual parameters.
5.	Select the parameter (1) capacity.
6.	Confirm by pressing the  button Enter capacity with the numeric keypad and subsequently confirm with  button.
7.	Use the  button to return to the parameter list
8.	Select the parameter (2) readability. Confirm by pressing the  button Enter readability via the numeric keypad and subsequently confirm with  button. (Possible values: 1,2,5,10,20,50)
9.	Use the  button to return to the parameter list

<p>10. Select the parameter (3) decimal point.</p> <p>Confirm by pressing the  button. Enter the decimal point (decimal places) via the numeric keyboard. Possible values: 0-3)</p> <p>11. Confirm by pressing the  button</p>
<p>12. Use the  button to return to the parameter list</p>
<p>13. Select the parameter (8) Mode. Enter the corresponding platform quantity via the numeric keyboard. (up to 6 platforms available)</p>
<p>14. Use the  button to return to the parameter list</p>
<p>15. Select the parameter (5) Calibration.</p>
<p>16. Adjusting the balance (see item 6.8)</p>

13 Declaration of Conformity

To view the current EC/EU Declaration of Conformity go to:

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