Scout[®] STX Series Balances

User Guide

EN

ES

FR

DE

IT

PT

HU

SV

DA

PL

CS

NL

1. SAFETY INFORMATION

Definition of Signal Warnings and Symbols

- WARNING** For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.
- CAUTION** For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.
- Attention** For important information about the product.
- Note** For useful information about the product.

Warning Symbols

General hazard

Electrical shock

Safety Precautions

CAUTION: Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the local AC power supply voltage is within the input voltage range printed on the AC adapter's ratings label.
- Only connect the AC adapter to a compatible grounded socket.
- Position the instrument such that the AC adapter can be easily disconnected from the socket.
- Position the power cord so that it does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in the user instructions.
- Do not operate the equipment in hazardous or explosive environments.
- Disconnect the equipment from mains power before cleaning or servicing.
- Service should only be performed by authorized personnel.

Intended Use

Use the instrument exclusively for weighing as described in the operating instructions. Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.

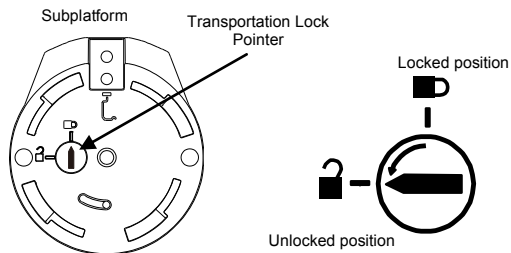
2. INSTALLATION

2.1 Selecting the location

The location must be sturdy, flat and level. Avoid locations with excessive air current, vibrations, heat sources or rapid temperature changes. Allow sufficient space around the instrument.


2.2 Releasing the Transportation Lock

On top of the balance, turn the pointer 90 degrees counter-clockwise to unlock.


2.3 Installing the Weighing Pan

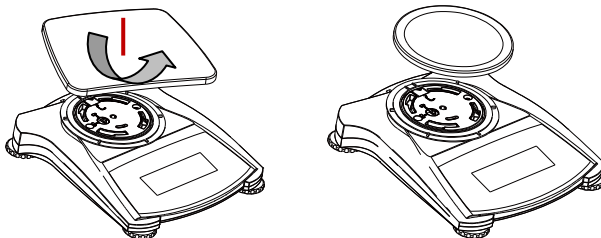Balances with a rectangular pan are placed into the sub-pan as shown and rotated counter-clockwise until it locks. Round pans are placed straight down on sub-pan.

2.4 Levelling the Instrument

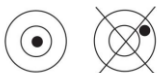To level the instrument, adjust the feet so the bubble is centered in the circle of the level indicator. Be sure the equipment is level each time its location is changed.



2.5 Connecting Power

AC Adapter Installation

Connect the AC adapter power cord to the instrument's power input connector, and then connect the AC plug to a suitable electrical outlet.

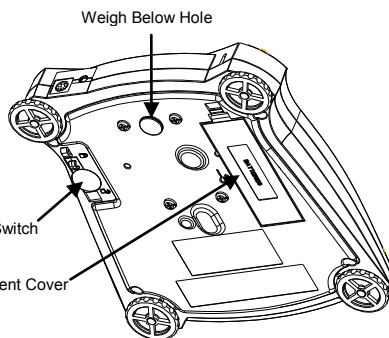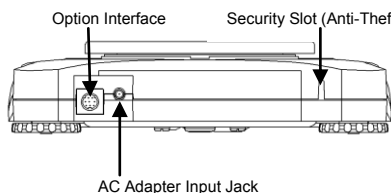Attention: Only use a AC adapter specified by OHAUS.

Battery Installation

Install the four "AA" batteries with polarity as shown in the battery compartment. When operating on batteries, long press the power button for up to 10 seconds to power on.

2.6 Initial Calibration

For best results, the instrument's calibration should be checked prior to first use. If adjustment is needed, refer to the Calibration section of the instruction manual.

3. OPERATION

3.1 Controls

Button	ZERO	TARE
Primary Function (Short Press)	On Turns the balance on Zero If balance is On, sets Zero	Tare Enter/clear a tare value.
Secondary Function (Long Press)	Off Turns the balance Off.	

3.2 Overview of Display, Home Screen

This equipment utilizes a touch-sensitive display. *Touch* areas and Buttons to control the equipment's functions.


Main Application Screen

Application		Capacity and readability
Instructional Messages	Place sample on the pan to begin averaging.	Result Field: Information varies by application Touch g to change unit
Stability (*), Net (>0<) indicators	* >0<	
Reference Fields	Gross: 0.000 g Averaging Time: 5 s Tare: 0.000 g Automatic Mode: On	Application Buttons: Function vary by application

3.3 MENU & SCREEN NAVIGATION

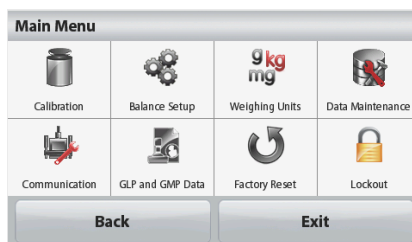Touch **Menu** to open the menu list.

-  **Calibration:**
Touch to view calibration options.
-  **Balance Setup:**
Touch to view balance settings.
-  **Weighing Units:**
Touch to view weighing units.
-  **Data Maintenance:**
Touch to view data maintenance options.
-  **Communication:**
Touch to view COM Device Settings and Print Settings.
-  **GLP and GMP Data:**
Insert user data for traceability.

-  **Factory Reset:**
Touch to do a Factory reset.
-  **Lockout:**
Touch to access the Lockout menu list.

3.3 Using the Balance

Weighing Application

1. If required, place an empty container on the pan and press **Tare**.
2. Add sample to the pan or container. The display shows the weight of the sample.

4. MAINTENANCE

4.1 Calibration

When the Balance is first installed, and when it is moved to another location, it must be calibrated to ensure accurate weighing results.

4.2 Cleaning

WARNING: Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.

Attention: Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The exterior surfaces of the instrument may be cleaned with a cloth dampened with water and a mild detergent.

4.3 Troubleshooting

For technical issues contact an Authorized Ohaus Service Agent. Please visit our website www.ohaus.com to locate the Ohaus office nearest you.

5. TECHNICAL DATA

The technical data is valid under the following conditions:

- Indoor use only
- Altitude: Up to 2000 m
- Operating temperature range: +5 °C to +40 °C
- Specified temperature range: +10 °C to +40 °C
- Relative humidity: 10% to 80% at 31°C, decreasing linearly to 50% at 40°C, non-condensing
- Mains supply voltage fluctuations: up to ±10% of the nominal voltage
- Installation category II
- Pollution degree: 2

SPECIFICATIONS:

MODEL	STX123	STX223	STX222	STX422	STX622	STX1202
Capacity	120 g	220 g	220 g	420 g	620 g	1000 g
Readability d	0.001 g	0.001 g	0.01 g	0.01 g	0.01 g	0.01 g
Repeatability (std. dev.)	2 d	2 d	1 d	1 d	1 d	2 d
Linearity	3 d	3 d	1 d	1 d	2 d	3 d
Power	AC power adaptor input 100-240 V 50/60 Hz and output 5V/1A DC, or 4 AA batteries					

MODEL	STX2202	STX421	STX621	STX2201	STX6201	STX8200
Capacity	2200 g	420 g	620 g	2200 g	6200 g	8200 g
Readability d	0.01 g	0.1 g	0.1 g	0.1 g	0.1 g	1 g
Repeatability (std. dev.)	2 d	1 d	1 d	1 d	1 d	1 d
Linearity (g)	3 d	1 d	1 d	1 d	2 d	1 d
Power	AC power adaptor input 100-240 V 50/60 Hz and output 5 V/1A DC, or 4 AA batteries					

6. COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	This product complies with the EU Directives 2014/30/EU (EMC), 2014/35/EU (LVD) and 2014/31/EU (NAWI). The EU Declaration of Conformity is available online at www.ohaus.com/ce .
	This product complies with the EU Directive 2002/96/EC (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to www.ohaus.com/weee .
	AS/NZS 61000.6.1, AS/NZS 61000.6.3

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P/N 30269065 C © 2021 Ohaus Corporation, All rights reserved
Ohaus Corporation